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1. Santos J, Mendiola JA, Oliveira MBPP, Ibanez E, Herrero M. Sequential determination of fat- and water-soluble vitamins in green leafy vegetables during storage. *J. Chromatogr. A*. 2012; 1261:179-188.
2. <http://www.prweb.com/releases/2011/10/prweb8906640.htm>
3. Blake CJ. Status of Methodology for the Determination of Fat-Soluble Vitamins in Foods, Dietary Supplements, and Vitamin Premixes, *J. AOAC Inter*. 2007; 90(4):897-910.
4. Gomis DB, Fernandez MP, Gutierrez Alvarez MD. Simultaneous determination of fat-soluble vitamins and provitamins in milk by microcolumn liquid chromatography, *J. Chromatogr. A*. 2000; 891:109-114.
5. Salo-Vaananen P, Ollilainen V, Mattila P, Lehtikoinen K, Salmela-Molsa E, Piironen V. Simultaneous HPLC analysis of fat-soluble vitamins in selected animal products after small-scale extraction, *Food Chem*. 2000;71:535-543.
6. Aubin A. Analysis of fat-soluble vitamin capsules using UltraPerformance Convergence Chromatography. Waters Application Note 720004394en. 2012 June.
7. Nelis HJCF, De Roose J, Vandenbaviere J, De Leenheer AP. Non-aqueous reversed-phase liquid chromatography and fluorimetry compared for determination of retinol in serum. *Clin. Chem*. 1983; 29(7):1431-1434.
8. Parriss NA. Non-aqueous reversed phase liquid chromatography: a neglected approach to the analysis of low-polarity samples. *J Chromatogr*. 1978; 157:161-170.

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©2013年沃特世公司 中国印制